



Note: Read the corresponding technical documentation for handling and safety reasons.


AZURA® inline check valve

Valid for order no. AZZ00FB & AZZ00FC
(spring-loaded)

Install 1/16" capillaries with a fitting (not included) with an UNF 10-32 thread to the inline check valve ① / ③. The arrow indicates the direction of the flow.



Note: The notch at the check valve ② indicates the inlet side.

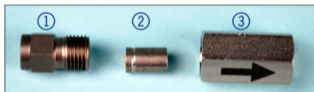

Exchanging the check valve

1. Unscrew inlet bushing ① from filter holder ③.
2. Exchange valve* ②.
3. Reattach the bushing ①.
Max. torque is 5 Nm.

* AZZ00FB: Order no. for spare valve is A06841
AZZ00FC: Order no. for spare valve is A068411